AGITADOR ORBITAL DIGITAL NAHITA BLUE

NAHITA BLUE DIGITAL ORBITAL SHAKER

Modelo / Model: 685/2

Referencia / Code: LJN001

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Key features

- Brushless motor for durability, stable, quiet and maintenance-free operation.
- LCD display for speed and time indication.
- The upper part of the motor is provided with a special waterproof insulation device, which completely solves the problem of liquid leakage and prevents the equipment from damage.
- The work tray is made of durable, corrosion-resistant engineering plastic. Four special screws are used to secure the front and rear of the tray, which facilitate disassembly and cleaning.

Specifications

- Display: LCD type, speed and time indication, memory of operation parameters (last used settings)
- Work tray: 300 x 300 mm, with anti-slip rubber and two elastic bands
- Rotation: Horizontal 360°
- Amplitude: 24 mm
- Speed range: 30-240 rpm
- Timing range: 99h 59 min 59 s
- Maximum shaking weight: 15 kg
- Motor: Brushless, maintenance-free operation
- Acoustic alarm at the end of shaking
- Power: 40W
- Ambient working temperature: -5 to 40°C
- Power supply: 220VAC, 50Hz
- Aprox. size: 370x300x150 mm
- Weight: 11 kg

Usage method

1. Speed setting: Turn the SPEED control knob to the required speed. Press the START/PAUSE key to start the agitation.
To finish press the STOP key.
During operation the SPEED key is locked, and it is not possible to change the speed. If you need to change the speed, first press SPEED to unlock and then adjust the speed.
2. Timer function: Press the TIME button to set the operation time.
This time is set in hours, minutes and seconds. When the number flickers, turn the TIME knob and set the desired value. Wait until the next number flickers to set its corresponding value.
You can also change from hours to minutes and seconds by pressing TIME.
If you do not want to set an operation time (continuous operation) set the time to 00:00:00.

3. Disassembly of the working tray: Loosen the four bakelite screws and remove the tray for cleaning.

Maintenance and cleaning

Proper use and maintenance will keep the equipment in good working condition and extend its useful life.

In routine work, keep the device clean and dry. If liquid spills on the working tray, wipe it up in time to prevent it from leaking into the agitator. If liquid enters the unit, the power supply must be shut off immediately.

Use a non-abrasive cleaning solution to clean the external surface. This solution should not seep into the interior of the shaker. Do not connect the power supply until the entire surface is dry.

Before maintenance or cleaning, disconnect the unit from the power supply.